

PRODUCT DESCRIPTION

PRODUCT NAME: GLYCOSEP™ C GUARD COLUMN
PRODUCT CODE: GKI-4721G
STORAGE: Room Temperature

Guard column specifically for use with GKI-4721 GlycoSep C HPLC Column (PPC 149). The GlycoSep C Guard Column is used in series with the GlycoSep C HPLC Column to remove sample impurities.

DIRECTIONS FOR USE:

Connect the inlet end of the GlycoSep C Guard Column (see figure below) to the HPLC system after the injector and before the GlycoSep C HPLC Column. Tighten the connection with fingers only; do not use a wrench.

Before connecting the GlycoSep C HPLC Column, run the mobile phase to purge air completely from the GlycoSep C Guard Column and the tubing which will connect to the GlycoSep C HPLC column.

Stop the pumps and connect the GlycoSep C HPLC Column downstream of the guard column. Make sure the flow indicators on both the guard column and the GlycoSep C HPLC Column are aligned in the proper direction.

Use the guard column as described in the product literature provided with the GlycoSep C HPLC Column.

NOTE: Minimize the tubing length between the injector and detector to reduce mixing effects.

