



## PRODUCT DESCRIPTION

PRODUCT NAME: GLYCOSEP™ R GUARD COLUMN  
PRODUCT CODE: GKI-4727G  
STORAGE: Room Temperature

Guard column specifically for use with GKI-4727 GlycoSep R HPLC Column (used in series to remove sample impurities).

### DIRECTIONS FOR USE:

Connect the inlet end of the guard column to the HPLC system after the injector and before the GlycoSep R column. Tighten the connection with fingers only; do not use a wrench.

With the R column disconnected, run the mobile phase to purge air in the guard column and the tubing to the R column.

Stop the pumps and connect the GlycoSep R column downstream of the guard column. Make sure the flow indicators on both the guard column and the R column are aligned in the proper direction.

Use the guard column as described in the product literature provided with the GlycoSep R column.

*NOTE: Minimize the tubing length between the injector and detector to reduce dispersion effects.*