



## PRODUCT DESCRIPTION

PRODUCT NAME: GLYCOSEP™ N GUARD COLUMN  
PRODUCT CODE: GKI-4728G  
STORAGE: Room Temperature

Guard column specifically for use with GKI-4728 GlycoSep N HPLC Column (used in series to remove sample impurities).

### DIRECTIONS FOR USE:

Connect the inlet end of the guard column to the HPLC system after the injector and before the GlycoSep N column.

With the N column disconnected, run the mobile phase to purge air in the guard column and the tubing to the N column.

Stop the pumps and connect the GlycoSep N column downstream of the guard column. Make sure the flow indicators on both the guard column and the N column are aligned in the proper direction.

Use the guard column as described in the product literature provided with the GlycoSep N column.

*NOTE: Minimize the tubing length between the injector and detector to reduce dispersion effects.*